

Fusion BioLabs X24 (Linear 24-amino acid peptide) Peptide Phage Display Library Information

SKU# PDL-04: Ready-to-Use X24 Linearized Peptide Phage Display Library Kit

Synthetic Format: Randomized 24-amino acid peptide with 6xHis Tag and HA Tag (OmpA leader sequence- **X24** peptide -6xHis tag-HA tag- Amber stop codon)

Diversity: 8.2×10^{10}


The diversity and in-frame of the library was checked by NGS.

The **X24** phagemid-based phage display Library is a linear randomized segment of 24-amino acid peptides.

Kit contents

The following components are included in the Kit.

Component	Quantity	Composition
X24 peptide phage display library (ready-to-panning; 8.2×10^{12} pfu/ml)	1.0 ml	1×PBS with 50% glycerol
M13KO7 Helper Phage (2×10^{12} pfu/ml)	0.5 ml	1×PBS with 50% glycerol
Chemically Competent TG1 E. coli	0.5 ml	2×YT with 25% Glycerol
M13 Reverse Primer (1.6 μ M)	0.2 ml	1×TE Buffer

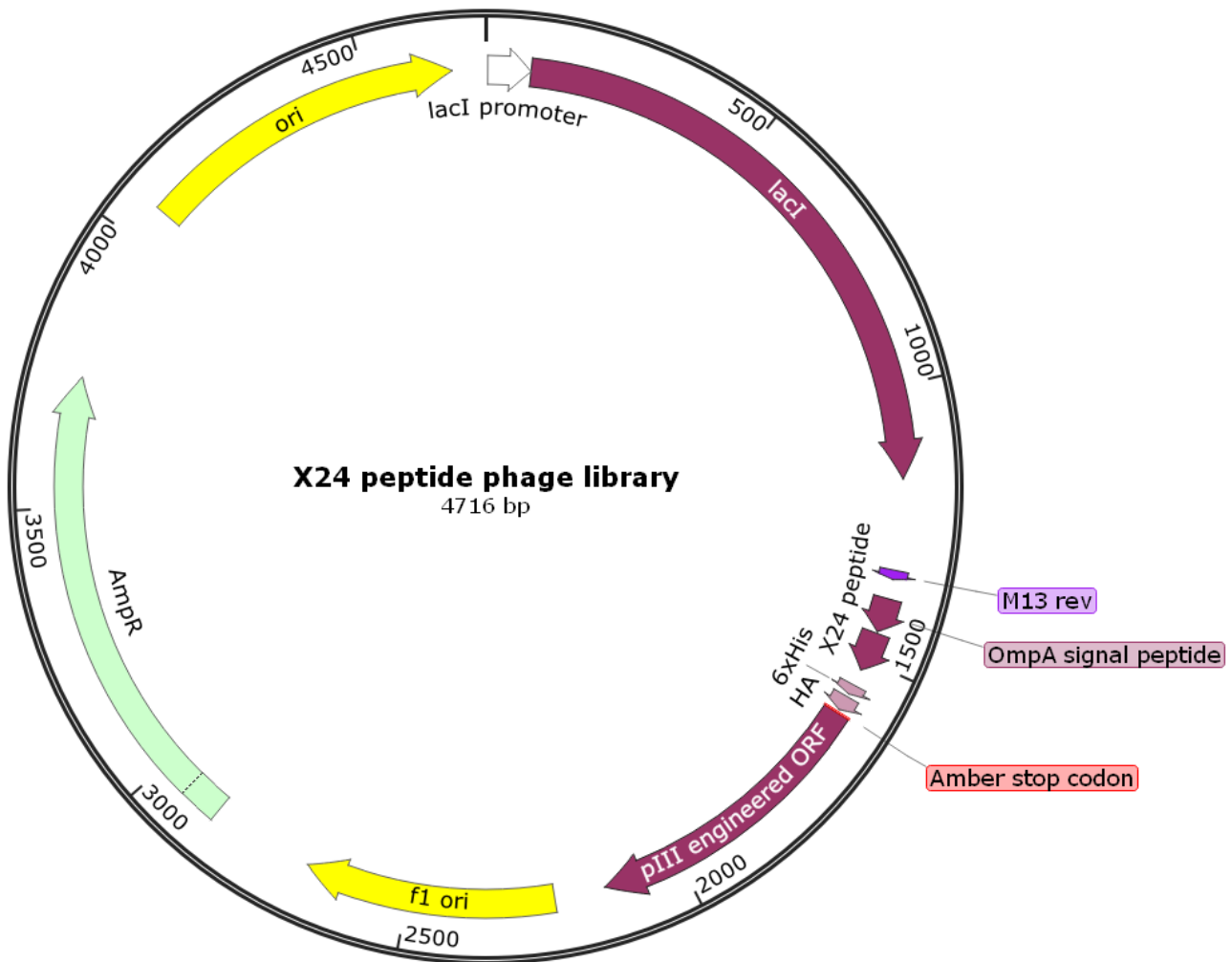

Comments for Fusion BioLabs Library Vector X24 Peptide Phage Display Library (one typical clone)

lac I promoter: bases 5-82

lac I ORF: bases 83-1165

M13 Reverse priming site: bases 1331-1347

OmpA leader sequence: bases 1383-1445

Polyhistidine (6xHis) site: bases 1551-1568

HA Hemagglutinin A (HA) epitope: bases 1575-1601

Amber stop codon: bases 1605-1607

pIII engineered ORF: 1608-2141

f1 origin: bases 2234-2689

Ampicillin resistance ORF: bases 2876-3736

pUC origin: bases 4068-4656